


## Centrifuge Tube (15 ml, 50 ml)

PakGent centrifuge tubes are made of high-purity polypropylene and molded using our advanced injection molding process to achieve efficient and stable centrifugation. They come with a recyclable plastic rack, which is environment friendly and clean. With a larger writing area, and special process to make our tubes excellent alcohol-resistant performance.




### Product Characteristics

Features	Description
Material	Polypropylene (PP), USP class IV
Centrifugal Force	Conical bottom: 15,000 xg, Conical bottom with skirt: 3,500 xg
Recommended Temperature Range	Autoclavable (121 °C) and freezable (-80 °C)
Sterility Level	SAL=10 <sup>-6</sup>
Cap Color	Dark blue, light blue caps available

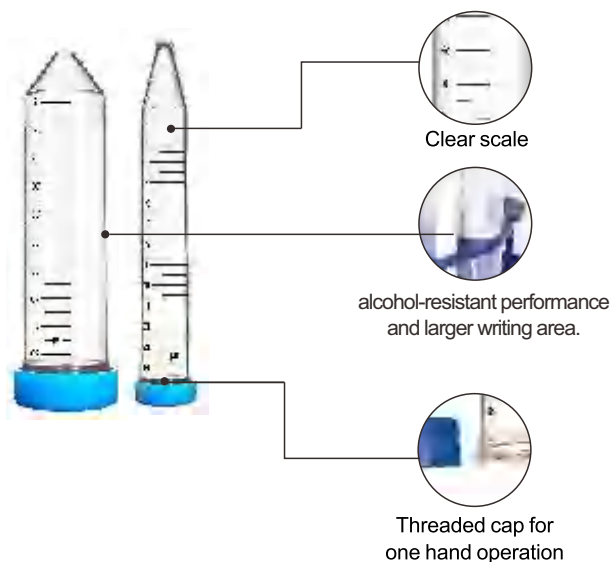

Uniform wall thickness and stable centrifugal effect.

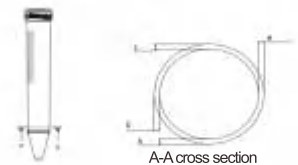

Zip bag design, more convenient for laboratory use.




Ergonomic, collapsible storage racks made from recycle material are better for lab and environment.

Available for both 15 mL and 50 mL sizes.



## Product Size Information

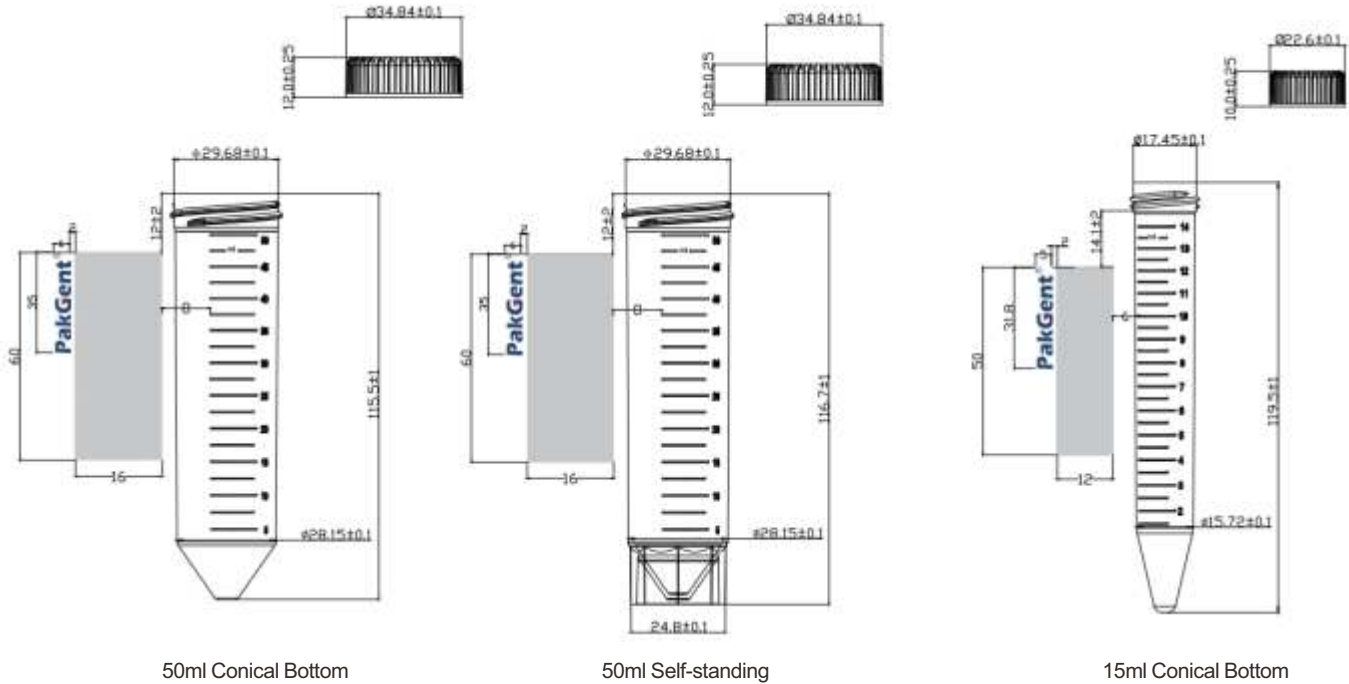

## Kindly Reminder

- After testing, the PakGent centrifuge tube graduation is accurate, but we still strongly recommend that the centrifuge tube graduation is only used as a reference . If the test requires accurate more precise volume measurement, it is recommended to use professional measuring tools.
- In order to ensure the centrifugal performance, it is recommended that the liquid should not exceed 80% of the max capacity when using a centrifuge tube for high-speed application.



It is recommended that the volume should not exceed 40 ml during centrifugation.



It is recommended that the volume should not exceed 12 ml during centrifugation.

- It is recommended to open the cap during the autoclave process to effectively prevent tube deformation of the tube.
- When storing at low temperatures, it is recommended to wrap sealing film around the connection between the tube and the cap to ensure better sealing function.

## Ordering information

Part No.	Description	Package
CT50R-L	50 ml conical bottom, alcohol-resistant performance, light blue cap, rack pack, 15,000 xg	25 pcs/rack, 20 racks/case
CT50B-L	50 ml conical bottom, alcohol-resistant performance, light blue cap, bulk pack, 15,000 xg	25 pcs/bag, 20 bags/case
CT50RS-L	50 ml conical bottom with skirt, alcohol-resistant performance, light blue cap, rack pack, 3,500 xg	25 pcs/rack, 20 racks/case
CT50BS-L	50 ml conical bottom with skirt, alcohol-resistant performance, light blue cap, bulk pack, 3,500 xg	25 pcs/bag, 20 bags/case
CT15R-L	15 ml conical bottom, alcohol-resistant performance, light blue cap, rack pack, 15,000 xg	25 pcs/rack, 20 racks/case
CT15B-L	15 ml conical bottom, alcohol-resistant performance, light blue cap, bulk pack, 15,000 xg	25 pcs/bag, 20 bags/case



# PakGent®

Save every drop of precious reagent for you.

**PakGent Bioscience (Suzhou) Co., Ltd.**  
Tel: +86 (0) 512 6670 0017  
Address: Floor 1-3, Building 33, Alloy Industrial Park, Wuzhong District, Suzhou, China.

**PakGent Bioscience (Changshu) Co., Ltd.**  
Tel: +86 (0) 512 5298 8330  
Address: No.9 Lecheng Road, Bixi Street, Changshu, Jiangsu Province, China.

**PakGent Bioscience (Shanghai) Co., Ltd.**  
Tel: +86 (0) 512 6670 0017  
Address: Building C, No 888, Lake west road, Lin gang section, Free Trade Zone, Shanghai, China.